


QUICK INFO



How to activate the Fridge-tag®



You receive the device in "sleep mode". The LCD screen is blank.

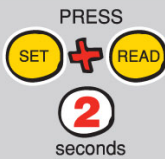

To activate the device press the **SET** and the **READ** buttons simultaneously for more than 2 seconds. Four zeros will appear on the screen. The first zero will flash.

Setting the time

Use the **READ** button to adjust the time. Each time you press the **READ** button, the number in the flashing digit will increase by 1. For example, if you want to set the time to 12:42 you have to perform the following steps:



Press **READ**. The first digit is flashing: Press **READ** once. "1" will appear as the first digit.



Press **READ**. The second digit will start flashing. Press **READ** two times, when "2" appears as the second digit.

Setting the time



Press **READ**. The third digit will start flashing. Press **READ** four times to set the digit as "4".



Press **READ**. The third digit will start flashing. Press **READ** two times to set the digit as "2".



Press **SET** to save. If during this operation, you press **READ** more then you were supposed to, continue pressing the **READ** button until you obtain the desired number, then press **SET** button to save your settings.

Setting the temperature measurement unit



When you finish setting the time, the "°C" sign will appear and blink at the right bottom corner. If you would like to record and read temperatures in **Celsius**, press **SET**.

If you want to record in **Fahrenheit**, press **READ** once and the "°F" sign will appear, then press **SET** to complete the activation.



The display now indicates the word **Loc** for a period of **10 minutes**, the **Fridge-tag®** does not measure the temperature to avoid false readings.

This allows an adaption to the changing environmental temperature when the device is placed in the refrigerator before normal recording starts. Place your device in the refrigerator now. The same lock function takes place after every reading (**by READ button pressing**) of the device.



Changing the time setting

If you made a mistake or wish to change the time setting, it can be readjusted. To adjust the time after the device is activated, you need to press and hold the **SET** button and then press the **READ** button shortly. Then you need to repeat the steps as described in "Setting the time". Additional time adjustments have no effect on the recording. Once the device is activated, it cannot be stopped. Adjustments can only be made for time setting and for changing the temperature measurement unit. After any correction, the adjustment mode is locked for **24 hours following** the midnight date shift. Thus, a new setting is possible only after the 2nd date shift again.

